

Storage Tanks and Accessories

Storage with a Difference



Guarantee the purity of your stored water

Pure water requires a storage system to prevent the degradation of your water quality. EMD Millipore's 30-, 60-, and 100-liter polyethylene (PE) storage tanks are designed to maintain consistent purity of stored water and provide effective protection against airborne contaminants.*

Prevent contamination

Water stagnancy can cause bacterial proliferation. Our optimal Automatic Sanitization Module (ASM) provides the ideal solution for the prevention of bacterial growth and biofilm formation on the inner surface of the storage tank. In addition, our advanced vent filter protects pure water from airborne contamination.

Distribute your stored water where it is needed

To provide pure water for use with all of their applications, laboratories need to be able to distribute stored water from their water purification system storage tanks.

- For distribution of **non-pressurized pure water**, a valve is conveniently located on the front of the EMD Millipore storage tanks.
- For convenient **distribution of pressurized pure water** from the storage tank, an E-POD® point-of-delivery dispenser can be connected to an Elix® Advantage or Milli-Q® Integral water purification system.
- For **automatic feed** of pure water, distribution valves on the base of the storage tank allow connection to other laboratory equipment such as glassware washing machines. Distribution pumps are also available if needed.

* A complete line of storage tanks is available, ranging in capacity from a few liters to several hundred liters. Your nearest EMD Millipore office will be able to guide you in the choice of the tank best suited to your needs.

Optimized pure water storage

The main concern when storing pure water is degradation of water purity over time. Only a strict choice of storage tank materials, associated with a careful design and appropriate protection against airborne contaminants, can ensure consistent water quality during storage.

Innovative storage tank design

EMD Millipore 30-, 60-, and 100-liter polyethylene storage tanks incorporate the latest technical developments and advanced features for stored water of consistent purity.

All tanks have a small footprint and are designed for wall-mounting if required. Underbench installation is also possible for some models.

Unique features

- Polyethylene selected for its minimum release of extractables
- Opaque walls block sunlight to prevent algae development
- Smooth inner surface prevents biofilm formation
- Cylindrical shape minimizes surface area in contact with water
- Conical bottom allows complete draining for cleaning and rinsing
- Pure water smoothly fed in at the bottom of the tank prevents absorption of carbon dioxide
- Front valve enables manual dispense of pure water
- Distribution valves permit connection to other laboratory equipment
- Hermetically sealed lid blocks air from entering the tank
- Large top opening allows manual cleaning during sanitization procedure
- Compact space-saving design


Fail-safe protection

- Sensor rod float switch system for automatic storage tank refill and indication of water level (% full)
- Overflow connected to the drain, in the unlikely event of a water system malfunction
- Direct display of stored water level on water purification system units
- Water sensor

Storage tanks designed for efficiency

Designed for efficiency

1. Hermetically sealed lid
2. Sensor rod float switch
3. Sanitary overflow
4. Blow-molded storage tank
5. Cylindrical shape
6. Conical bottom with distribution valves
7. Front dispensing valve
8. Advanced vent filter

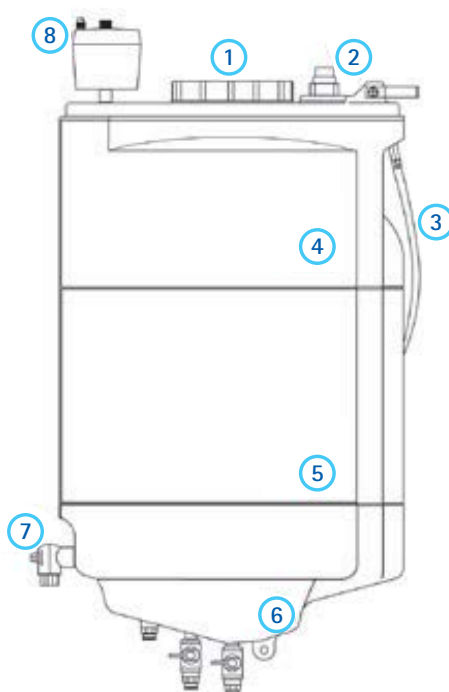

For details of the tests performed during the storage tank development process, please request the publication "R&D Notebook 1: Optimizing the storage of purified water for laboratory applications" (Ref. No.: RD001EN00) from your local EMD Millipore representative.

Storage Tank Accessories

In order to help ensure optimum purity and distribution of your stored water, EMD Millipore offers a range of accessories and connections for your storage tank, including the following items:


- Advanced Vent Filter
- Automatic Sanitization Module
- Air Gap Device
- E-POD® pure water remote dispenser
- Washer Distribution Kit
- Water Sensor
- Lab Close Kit

Advanced vent filter: important protection against airborne contaminants

Air is contaminated by carbon dioxide, particles, microorganisms, and volatile organic compounds that come mainly from the laboratory atmosphere. To protect pure water from all these contaminants, EMD Millipore has developed an advanced storage tank vent filter that includes:

- Activated carbon to adsorb volatile organics (including lab solvents such as acetone, chloroform, and methanol)
- A soda-lime bed to remove CO₂
- A Durapore® hydrophobic membrane for particle and bacteria retention

This advanced vent filter is recommended for the protection of high-resistivity water, such as Elix® product water, during storage. To protect RiOs™ reverse osmosis-quality water, a Durapore® 0.45 µm hydrophobic membrane vent filter is also available.

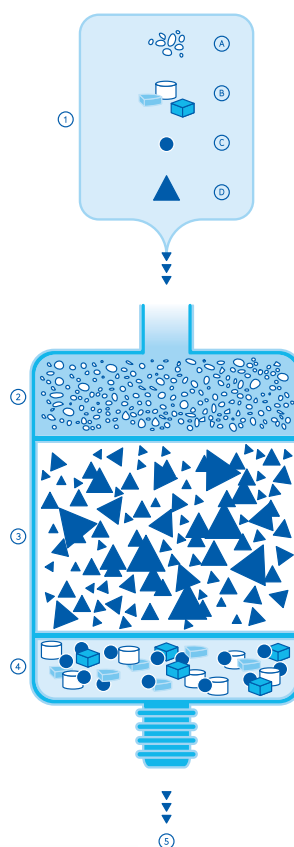

Advanced storage tank vent filter


1. Airborne Contaminants
 - A. Volatile Organics
 - B. Particles
 - C. Bacteria
 - D. CO₂
2. Volatile Organics absorption
3. CO₂ removal
4. Particle and Bacterial retention
5. Storage tank inlet
Purified air enters the storage tank



Automatic Sanitization Module (ASM): say "No!" to bacterial proliferation


Maintaining high purity water with low bacteria levels during storage is critical. If left to proliferate, trace levels of microorganisms present in pure water compromise water purity. This bacterial contamination is responsible for the formation of a biofilm — an accumulation of organic material made up of active and dead organisms, on the inner walls of the storage tank.

Even though chemical sanitization and mechanical scrubbing may be periodically performed, this biofilm is difficult to remove and is a frequent source of recontamination in stored water.

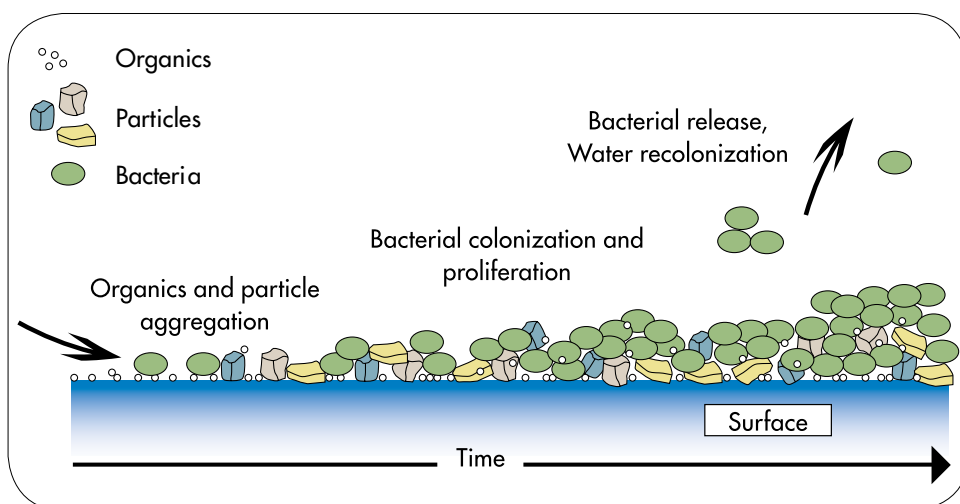

Biofilm formation over time

EMD Millipore's ASM is designed to prevent the growth and proliferation of bacteria and the resulting biofilm on the inner surface of EMD Millipore PE storage tanks. The ASM makes use of the germicidal properties of an ultraviolet (UV) light at 254 nm, which is fitted inside the tank.

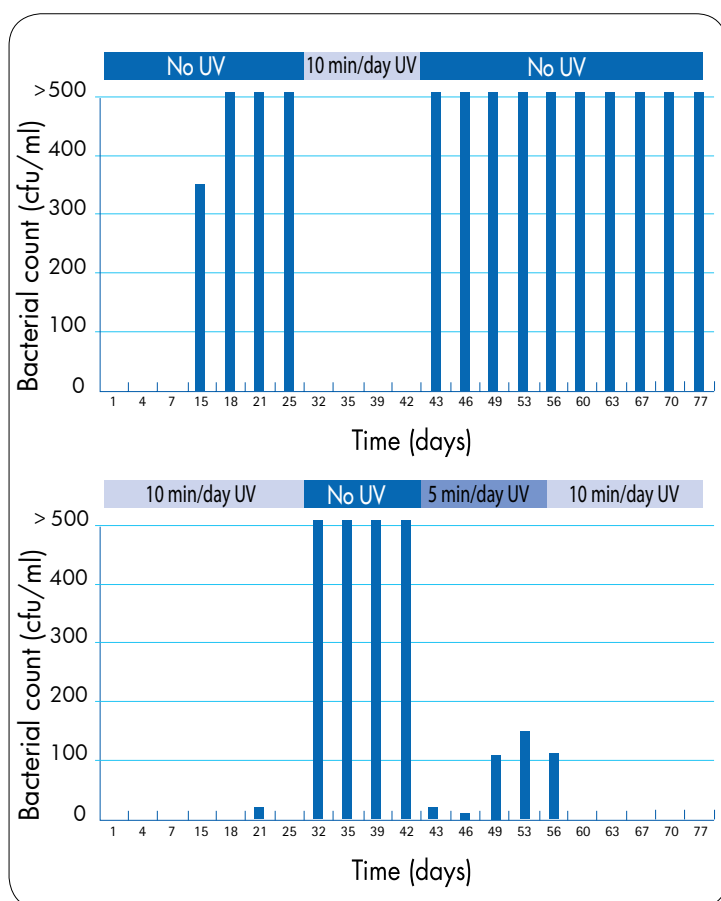

Germicidal effectiveness of the UV lamp

The ASM provides full flexibility for guaranteed results

- 254 nm UV lamp; selected for its germicidal effectiveness
- Pre-programmed intervals of 10-min / day automatic UV illumination for optimized efficiency
- Additional programmable and manual UV exposure possible to meet critical application requirements
- Up to 45 min /day of UV exposure for total flexibility
- Program daily time settings, UV cycles, and UV lamp operation displayed on the Millitrack® e-Solution dashboard
- UV lamp exchange alarm for easy maintenance
- Compact design allowing installation on top of the storage tank

10 minutes of daily UV exposure is sufficient

During development of the ASM, the UV lamp exposure cycles were optimized by examining the resulting bacterial reduction after exposure.

Two 60-liter storage tanks were fed by an intentionally contaminated reverse osmosis water purification system. The tanks were then emptied and refilled each day and alternately equipped with an ASM into which variable illumination times were programmed.

As shown in the graphs, 10 minutes per day of UV exposure were enough to make the tanks return to their original low bacterial levels.

Air gap device for protection against bacteriological contamination

Water purification systems and storage tanks sometimes require a connection to the drain.

Drains are typically dirty environments contaminated by microorganisms, and in particular, bacteria. Therefore, when the outlet of the water system reject tubing is pushed into the drain, there is a risk that bacteria could contaminate the inside of the reject tubing, and then progressively move to the water system.

One way to prevent this from occurring is to install an air gap device on the reject tubing. This allows the reject water flow to move through the tubing without touching the inside of the contaminated drain environment. Installing an air gap device is an easy and safe way to prevent the development of bacteria above the air gap level.



Air Gap Device

E-POD® pure water remote dispenser: pure water where you need it

The E-POD® Elix® water point-of-delivery unit can be connected to an Elix® Advantage pure water system or Milli-Q® Integral pure and ultrapure water system to dispense pure water wherever it is needed in the lab.

Advantages of the E-POD® remote dispenser include:

- Improved bacterial water quality (less than 0.1 cfu/mL, with final filter)
- Versatility enabling use for multiple applications or users when a Millipak® or Biopak® polisher is fitted to the dispenser outlet
- Volumetric dispensing to save time
- Ergonomic design and ease of use
- Flexibility, with installation of up to three E-POD® units per system
- Information at a glance thanks to the color backlit screen on the dispenser base
- Space-saving small footprint



E-POD® pure water remote dispenser




Distribution pumps to meet increased demands

Washer Distribution Kit

A key use of pure water is as feed to laboratory appliances such as glassware washers, autoclaves, sterilizers and weathering devices.

The EMD Millipore Washer Distribution Kit provides cost-effective and convenient distribution of pure water to common laboratory appliances, with flow rates between 12.5 – 13.5 L/min (at 2 bar or 30 psi, depending on voltage).

Installation of the small-footprint kit is fast, easy, and universal, with bench, underbench, or wall-mounted options. Users profit from a silent, automatic supply of pure water when required.

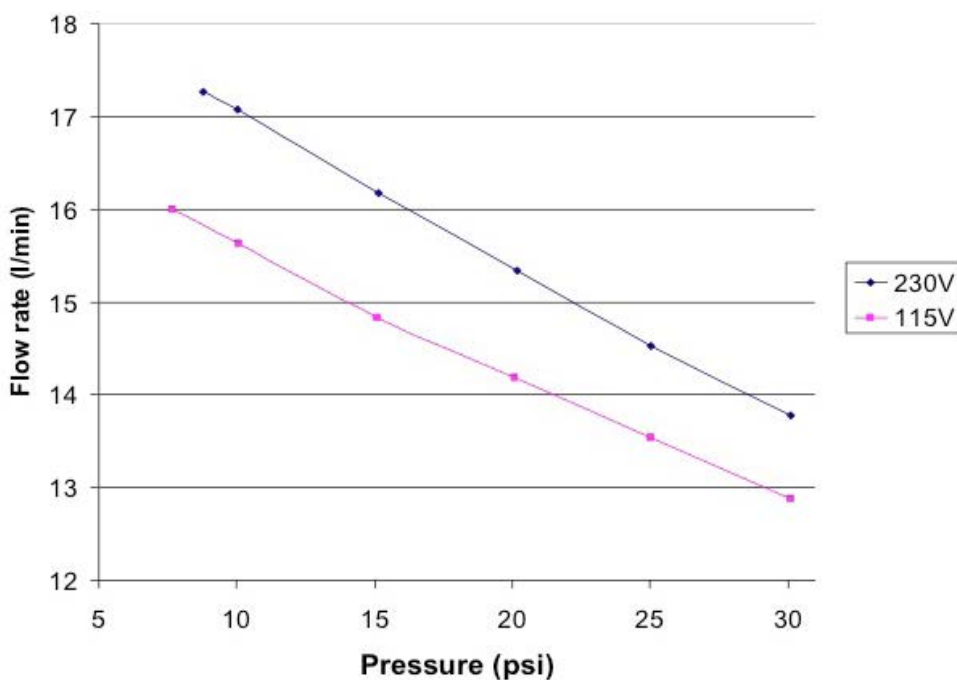

The graph shows characteristics of EMD Millipore distribution pumps. Pumps provide a long service lifetime and quiet operation.



Washer Distribution Kit

Water sensor for control over feed water supply

If there is water on the floor, the water sensor enables shutdown of the feed water supply in order to prevent a lab flood.

Lab Close Kit keeps your system in top condition when you're away

When your facility is closed for an extended time – such as vacation periods – the Lab Close Kit will avoid water purification system standstill during these long periods. The Lab Close Kit intelligently optimizes the consumption of water and electricity by your system, preventing the negative effects of nonuse, such as bacteria buildup. Your water purification system remains in top condition, ready for immediate use upon your return.

Specifications & Ordering Information

Specifications for Polyethylene Storage Tanks

There are several different storage tank sizes available in this range:

30-liter Storage Tank*

| | |
|-------------------------|--------------------|
| Diameter | 380 mm (14.82 in.) |
| Height | 600 mm (23.4 in.) |
| Maximum Usable Capacity | 25 L |
| Weight (full) | 30 kg (66.14 lb) |

60-liter Storage Tank

| | |
|-------------------------|--------------------|
| Diameter | 380 mm (14.82 in.) |
| Height | 840 mm (32.76 in.) |
| Maximum Usable Capacity | 54 L |
| Weight (full) | 59 kg (130.07 lb) |

100-liter Storage Tank*

| | |
|-------------------------|---------------------|
| Diameter | 380 mm (14.82 in.) |
| Height | 1255 mm (48.95 in.) |
| Maximum Usable Capacity | 91 L |
| Weight (full) | 98.5 kg (217.15 lb) |

* For 30-liter and 100-liter storage tanks, underbench models are also available.

Ordering Information

| Description | Catalogue No. |
|-------------|---------------|
|-------------|---------------|

Polyethylene Storage Tanks

| | |
|------------------------------|-----------|
| 30-liter PE tank | TANKPE030 |
| 30-liter PE underbench tank | TANKBI030 |
| 60-liter PE tank | TANKPE060 |
| 100-liter PE tank | TANKPE100 |
| 100-liter PE underbench tank | ZBITANK01 |

| Description | Catalogue No. |
|-------------|---------------|
|-------------|---------------|

Accessories

Advanced Vent Filter

| | |
|--|-----------|
| Advanced vent filter (for Elix® water purification systems) | TANKMPK01 |
| Standard vent filter (for RiOs™ water purification systems) | TANKMPK02 |

Automatic Sanitization Module

| | |
|---|-----------|
| Milli-Q® Integral; Milli-Q® Direct, Elix® Advantage, and Elix® Reference water purification systems | TANKASMIN |
| ASM for RiOs™ / Elix® / AFS® Essential water purification systems | TANKASMES |

Air Gap Device

| | |
|----------------|-----------|
| Air Gap Device | AIRGAP001 |
|----------------|-----------|

E-POD® Pure Water Remote Dispenser

| | |
|-----------------------------|-----------|
| E-POD® pure water dispenser | ZRXSP0D01 |
|-----------------------------|-----------|

Washer Distribution Kit

| | |
|---------------------------------------|-----------|
| Washer Distribution Kit (Left) 230 V | ZWDK5L100 |
| Washer Distribution Kit (Left) 115 V | ZWDK6L100 |
| Washer Distribution Kit (Right) 230 V | ZWDK5R100 |
| Washer Distribution Kit (Right) 115 V | ZWDK6R100 |

Water Sensor

| | |
|--------------------------------------|--|
| Connection from the system | ZFWATDET4 |
| Connection from the tap water source | ZFWATDET1 (120 V) or ZFWATDET2 (230 V) |
| Water sensor with cable | TANKLK002 |

Lab Close Kit

| | |
|---|-----------------------------------|
| Milli-Q® Integral; Elix® Advantage; Elix® Reference; and Milli-Q® Direct water purification systems | LABCLOSE1 |
| RiOs™ / Elix® / AFS® Essential water purification systems | Included, setting to be activated |

For more information, please visit our website:
www.emdmillipore.com/labwater

PB5506ENUS Copyright 2015 EMD Millipore Corporation, Billerica, MA., U.S.A.

EMD Millipore, the M mark, Milli-Q, Elix, AFS, E-POD, Millitrac, Millipak, Biopak, and Durapore are registered trademarks of, and RiOs is a trademark of, Merck KGaA, Darmstadt, Germany. All rights reserved.