


## PRECISION Standard Electronic Balances

MODELS TS120S, TS400S, TS400D, TS4KS and TS4KD

Instruction Manual

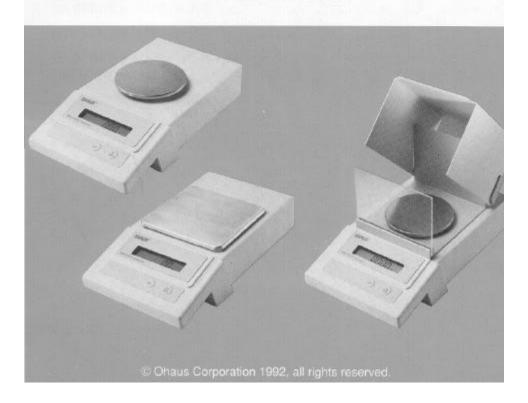

WARNING: THIS EQUIPMENT GENERATES, USES, AND CAN RADIATE RADIO FREQUENCY ENERGY AND IF NOT INSTALLED AND USED IN ACCORDANCE WITH THE INSTRUCTION MANUAL, MAY CAUSE INTERFERENCE TO RADIO COMMUNICATIONS. IT HAS BEEN TESTED AND FOUND TO COMPLY WITH CLASS "A" REQUIREMENTS IN BOTH PART 15 OF FCC RULES AND THE RADIO INTERFERENCE REGULATION OF THE CANADIAN DOC. THIS EQUIPMENT DOES NOT EXCEED THE LIMITS FOR RADIO NOISE EMISSIONS. OPERATION OF THIS EQUIPMENT IN A RESIDENTIAL AREA MAY CAUSE UNACCEPTABLE INTERFERENCE TO RADIO AND TV RECEPTION REQUIRING THE OPERATOR TO TAKE WHATEVER STEPS ARE NECESSARY TO CORRECT THE INTERFERENCE.

LE PRÉSENT APPAREIL NUMÉRIQUE N'EMET PAS DE BRUITS RADIOÉLECTRIQUES DEPASSANT LES LIMITES APPLICABLES AUX APPAREILS NUMÉRIQUES DE CLASSE A PRESCRITES DANS LE REGLEMENT SUR LE BROUILLAGE RADIOÉLECTRIQUE EDICTÉ PAR LE MINISTERE DES COMMUNICATIONS DU CANADA.

#### **TABLE OF CONTENTS**

| UNPACKING                | 4  |
|--------------------------|----|
| INSTALLATION             | 5  |
| Environment              | 5  |
| Draft Shield             | 5  |
| Pan and Pan Support      | 6  |
| AC Adapter               | 6  |
| Turning the Balance ON   | 7  |
| CHECKING CALIBRATION     | 7  |
| OPERATION                | 8  |
| Auto Range Models        | 8  |
| Averaging Level          | 8  |
| Weighing                 | 9  |
| Taring                   | 9  |
| RS-232 INTERFACE         | 10 |
| CALIBRATION PROCEDURE    | 10 |
| INTERNAL SWITCH SETTINGS | 13 |
| TROUBLESHOOTING          | 15 |
| SPECIFICATIONS           | 16 |
| CARE AND MAINTENANCE     | 16 |
| REPLACEMENT PARTS        | 17 |
| ACCESSORIES              | 17 |
| LIMITED WARRANTY         | 18 |

#### **PREFACE**


Your OHAUS® PRECISION *Standard* balance is a precision weighing instrument that is designed to be accurate and easy to operate. This manual explains how to use your balance properly and should be read before operation.

#### UNPACKING

Your PRECISION Standard balance was shipped with the following items:

- a pan
- · a pan support
- · an AC power adapter
- a draft shield kit (TS120S and TS400D only) includes draft shield and snap clamp
- · this instruction manual
- · your warranty card

It is recommended to save the carton and packing material for storing and/or transporting the balance.

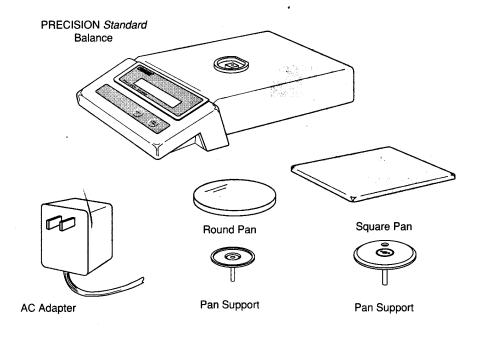

(Round pan and support for Models TS120S, TS400S, and TS400D)

(Square pan and support for Models TS4KS and TS4KD)

#### INSTALLATION

#### **Environment**

The balance should always be used in an environment which is free from excessive air currents, corrosives, vibration, and temperature or humidity extremes. These factors will affect displayed weight readings.

#### DO NOT install the balance:

- next to open windows or doors causing drafts or rapid temperature changes.
- · near air conditioning or heat vents.
- near vibrating, rotating or reciprocating equipment.
- near magnetic fields or equipment that generates magnetic fields.
- · on an unlevel work surface.

#### Draft Shield (TS120S and TS400D)


To install the draft shield:

 Position the draft shield on top of the balance as shown.
 Make sure the rib at the front of the draft shield base butts up against the raised lip of the display panel.

The snap clamp should be oriented so that the double clip is toward the front of the

balance.

Insert the double clip of the snap clamp into the opening in the draft shield base, and clip the draft shield base to the balance. Press the rear clip of the snap clamp into place.

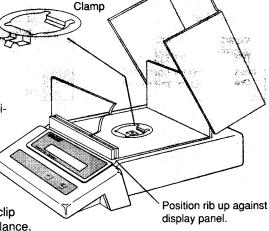


Snap

#### Pan and Pan Support

#### Square Pan

Insert the pan support into the hole in the weighing machanism as shown in the illustration. Make sure the hole in the pan support faces the rear of the balance. Once installed, the pan support should not rotate.


The pan has a guide pin which protrudes from the bottom. Place the pan on the support making sure the guide pin is inserted in the hole in the pan support.

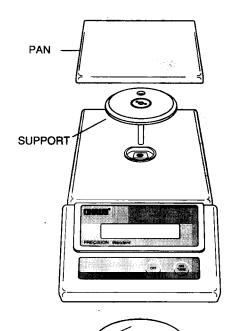

#### Round Pan


Insert the pan support into the hole in the weighing mechanism.

Place the pan on the support.

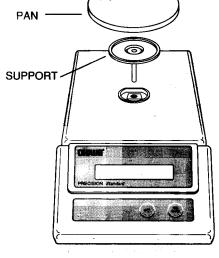

#### **AC Adapter**

Plug the molded connector of the adapter into the receptacle at the rear of the balance (see diagram). Plug the adapter into a convenient AC outlet.

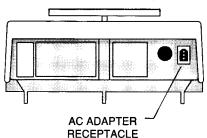

#### **Turning the Balance ON**

With no load on the pan, switch the balance ON by pressing the TARE button. When first switched ON, all segments of the display should be on as shown in the illustration.

This "display check" will be displayed briefly, then the model number of the balance will be displayed. If the balance is a dual range model (TS400D or TS4000D), the word "dUAL" will then be displayed. The display will momentarily blank and then indicate zero.

**\*.8.8.8.8.8** 9

£5400

#### CHECKING CALIBRATION

Before using the balance, calibration should be checked. The balance has been calibrated before shipment, however, it could be influenced by factors such as:

- variations in the earths gravitational field at different latitudes of the world
- rough handling
- · changes in work location


Weights required for checking calibration are listed in the adjacent table.

To check the balance's calibration, place the appropriate weight on the center of the pan and read the displayed weight.

If the displayed weight differs from the known weight by more than acceptable limits (see SPECIFICATIONS on page 16), the balance must be calibrated as explained in the "CALIBRATION PROCEDURE" section .

| CALIBRATION CHECK WEIGHTS |        |      |
|---------------------------|--------|------|
| MODEL                     | WEIGHT |      |
| TS120S                    | 100g   | . 75 |
| TS400S                    | 400g   | 14 · |
| TS400D                    | 400g   |      |
| TS4KS                     | 4kg    |      |
| TS4KD                     | 4kg    |      |

Weights must meet ASTM Class 1 Tolerance. Calibration weights are available as accessories (see page 17).

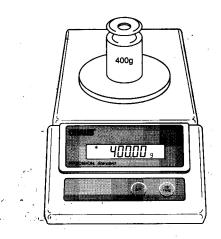

#### **OPERATION**

### Auto Range Models (TS400D and TS4KD)

Auto range balances offer both a fine range (lower capacity/higher readability) and a coarse range (higher capacity/lower readability). When first turned on, the balance is in the fine range. It remains in this range until the weight on the pan exceeds the fine range capacity. When weight on the pan is greater than the fine range capacity, the balance switches to the coarse range.

If weight on the pan falls back to within the fine range capacity, coarse range readability remains in effect until you tare the balance with no weight on the pan.

#### **Averaging Level**

Averaging level compensates for vibration or excessive air currents on the pan. During operation, the balance continually takes weight readings from the load cell. Successive readings are then digitally processed to achieve a stabilized display.

PRECISION Standard balances allow you to select a High or Low averaging level depending on how much processing you need to obtain a stabilized display:

HIGH more processing, greater stability, slower stabilization time.

LOW less processing, less stability, faster stabilization time.

**NOTE:** Averaging level does not affect balance accuracy.

The balance is shipped from the factory with Averaging Level set to Low. To change the averaging level, refer to the section titled "INTERNAL SWITCH SETTINGS".

Weighing

1. Press ON to rezero the display.

- 2. Place the object(s) or material to be weighed on the pan.
- 3. Wait for the stability indicator to appear before reading the weight.


# STABILITY INDICATOR \* 12345 9

#### **Taring**

When weighing material or objects that must be held in a container, taring enables you to store the container weight in the balance's memory, separate from the weight of the material in the container.


- 1. Place an empty container on the pan. Its weight will be displayed.
- 2. Press ON TARE

The display will show zero and the container's weight will be stored in memory.

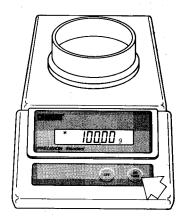

- Add material to the container. As material is added, its net weight will be displayed.
- Removing the container and material from the pan will cause the balance to display the container's weight as a negative number.

Tared weight remains in balance memory until ON is pressed again.

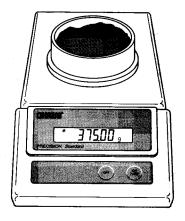

#### **RS-232 INTERFACE**

If your balance is equipped with the RS-232 Interface accessory, refer to the instructions provided with that accessory for operating information.

#### CALIBRATION PROCEDURE

PRECISION Standard balances can be calibrated in two ways: Span calibration or Linearity calibration.

**Span calibration** resets the balance's weighing range using two weight values: zero and a weight value at or near the balance's capacity.

**Linearity calibration** minimizes deviation between actual and displayed weights within the balance's weighing range. Three weight values are used: zero, a weight value within the balances weighing range, and a weight value at or near the balance's specified capacity.

Calibration should be performed as necessary to ensure accurate weighing. Weights required to perform the procedures are listed in the adjacent table.

NOTE: If calibration has been locked out, you will not be able to access it. Refer to the section titled "INTERNAL SWITCH SETTINGS" to enable it.

| CALIBRATION WEIGHTS |                      |                     |  |
|---------------------|----------------------|---------------------|--|
| MODEL               | LINEARITY<br>WEIGHTS | SPAN ONLY<br>WEIGHT |  |
| TS120S              | 50g,100g             | 100g                |  |
| TS400S              | 200g, 400g           | 400g                |  |
| TS400D              | 40g, 400g            | 400g                |  |
| TS4KS               | 2kg, 4kg             | 4kg                 |  |
| TS4KD               | 400g, 4kg            | 4kg                 |  |

Weights must meet ASTM Class 1 Tolerance. Calibration weights are available as accessories (see page 17).

### BEFORE BEGINNING CALIBRATION, MAKE SURE WEIGHTS ARE ON HAND.

If you begin calibration and realize weights are not available, either turn the balance off, or go through the procedure without weights. The balance will use previously stored calibration data

To start **SPAN** calibration, the BALANCE MUST BE **ON**.

Press and hold until "CAL" is displayed, then release it.

To start **LINEARITY** calibration, the BAL-ANCE MUST BE **OFF**.

Press and hold (N) until the balance turns on and "CAL" is displayed, then release it.

- When take is released "C 0 g" will be displayed indicating that no weight should be on the pan.
- With no weight on the pan,
   press ON TARE.


- The display will show "C" followed by the value of the weight which must be placed on the pan.
- Place the required weight on the pan and press ON TARE.


If Span calibration is being performed, proceed to step 7.

If Linearity calibration is being performed, continue with step 5.

[AL



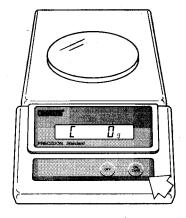

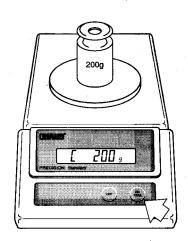

5. The display will show "- C -" momentarily, then "C" followed by the next weight to be placed on the pan.


**DO NOT disturb the balance** when "- C -" is displayed. Disturbances will result in improper calibration.

- 6. Place the required weight on the pan, then press (NARE).
- The display will show "- C -" while the balance recalibrates. When the weight on the pan is displayed along with the current unit indicator, the balance is recalibrated.

DO NOT REMOVE CALIBRATION WEIGHT(S) UNTIL WEIGHT VALUE IS DISPLAYED AND UNIT INDICATOR APPEARS.

#### **Calibration Protection**

Calibration may be "locked-out" to prevent unauthorized personnel from changing calibration. To lock out calibration refer to the section titled "DIP SWITCH SETTINGS".

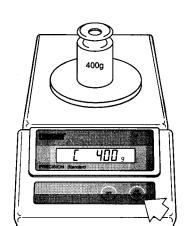

#### INTERNAL SWITCH SETTINGS

PRECISION *Standard* balances contain internal switches which may be set to do the following:


- Select a High or Low Averaging Level
- Disable/Enable calibration
- Disable/Enable the Print menu (only on balances equipped with the RS-232 option).


To access the switches, use the following procedure:

- 1. Disconnect power to the balance.
- 2. Remove the pan and pan support.
- 3. Using a philips screwdriver, remove the three screws which secure the cover to the balance. The balance will have to be turned over or on its side to access the screws as they are located on the bottom; two under the front panel and one on the rear (see diagram). HOLD THE COVER AND BASE TOGETHER WHILE REMOV-ING SCREWS. DO NOT ALLOW COVER OR BASE TO FALL. Remove the cover.

- 5. Locate the bank of four switches on the main circuit board under the top right portion of the display.
- 6. Set the desired switches using the following table as a guide.

| Function           | Switch | Setting                             |
|--------------------|--------|-------------------------------------|
| Averaging<br>Level | 1      | Closed = High<br>Open = Low         |
| Calibration        | 3      | Closed = Disabled<br>Open = Enabled |
| Print Menu*        | 4      | Closed = Disabled<br>Open = Enabled |

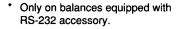

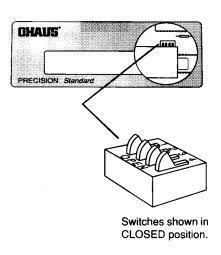

- 7. Replace the cover and fasten the three screws that were removed.
- 8. Replace the pan and pan support.
- 9. Reconnect power to the balance.

#### TROUBLESHOOTING

| SYMPTOM  | PROBABLE CAUSE(S)  | REMEDY  |
|--|--|---|
| Unit will not turn on.                         | Power adapter not plugged in or properly connected to balance. | Check power adapter connections.  |
| Incorrect weight reading.                      | Balance was not re-zeroed before weighing.                     | Press TARE with no weight on the pan, then weigh item.                              |
|  | Balance not properly calibrated.                               | Recalibrate correctly.  |
| Random segments displayed or display locks up. | Microprocessor lock-up.  | Turn balance off, then turn on again. If condition persists, unit must be serviced. |

#### **Error Codes**

- 2.0 Balance is unable to stabilize within time limit after taring. Environment is too hostile or balance needs recalibration.
- 3.0 Incorrect or no calibration weight used for calibration. Recalibrate with correct weights.
- 8.0 Hardware error causing an internal weight signal which is too low. Check if pan or pan support is off. If not, balance must be serviced.
- 8.1 Hardware error causing an internal weight signal which is too high. Have balance serviced.
- 8.2 Power-on load out of specification: Balance was turned on with load on pan or pan off balance. No load may be on pan when turned on and pan must be in place.
- 8.3 Overrange error. Load on pan exceeds capacity of balance.
- 9.8 Hardware error causing invalid calibration data checksum. Balance may need recalibration - particularly linearity calibration. If error persists, balance must be serviced.
- 9.9 Invalid temperature compensation checksum. Balance will work with default temperature compensation data. Error will re-occur each time balance is turned on. Have balance serviced.

#### **SPECIFICATIONS**

| MODEL                | TS120S                | TS400S        | TS400D             | TS4KS        | TS4KD    |
|----------------------|-----------------------|---------------|--------------------|--------------|----------|
| Capacity (g)         | 120                   | 400           | 400/40             | 4000         | 4000/400 |
| Readability (g)      | .001                  | .01           | .01/.001           | .1           | .1/.01   |
| Weighing Modes       |                       |               | grams only         |              |          |
| Tare                 |                       | Full Car      | pacity by Subtract | tion         |          |
| Precision/           |                       |               |                    |              |          |
| Reproducibility (g)  | .001                  | .007          | .007/.001          | .07          | .07/.01  |
| Linearity (g)        | .001                  | .01           | .01/.001           | .1           | .1/.01   |
| Sensitivity Drift    |                       |               |                    |              |          |
| (10 - 30 °C)         |                       |               | 10ppm/ °C          |              |          |
| Display              |                       |               | .6" LCD            |              |          |
| Stabilization Time   | 1 - 3 3 3 3 3 3       |               | 2.5 (sec)          |              |          |
| Operating            |                       |               |                    |              |          |
| Temperature          |                       | 50 -          | 104°F/10 - 40°C    | _            |          |
| Power Supply         | A                     | Adapter - 100 | , 120, 220, 240 V  | AC, 50/60 Hz |          |
| Calibration          | 1 264                 |               | Pushbutton         |              |          |
| Pan Size (in)        | 4.75 dia. 6.5W x 6.0D |               |                    |              |          |
| (mm)                 | 121 dia. 165W x 152D  |               |                    |              |          |
| Housing Size (in)    | 8.5W x 14.0D x 3.75H  |               |                    |              |          |
| (mm)                 |                       | 216           | SW x 356D x 95H    | -            |          |
| Net Weight (lb)/(kg) | , yes                 |               | 9.8/4.4            |              |          |

#### CARE AND MAINTENANCE

To keep the balance operating properly, the housing and pan should be kept clean and free from foreign material. If necessary, a cloth dampened with a mild detergent may be used. Keep calibration weights in a safe dry place.

#### SERVICE INFORMATION

If the Troubleshooting section does not resolve or describe your problem, you will need to contact an authorized Ohaus Service Agent. For Service assistance in the United States, please call Ohaus Corporation toll-free at (800) 526-0659. An Ohaus Product Service Specialist will be available to help you.

#### **REPLACEMENT PARTS**

| AC Adapters:                                | OHAUS<br><u>Part No.</u> |
|---|--------------------------|
| 100V  | 90524-11                 |
| 120V  | 90524-10                 |
| 220V  | 90524-13                 |
| 240V  | 90524-14                 |
| Pan - 4.75" dia.                            | 77262-10                 |
| Pan - 6.0" x 6.5"                           | 77298-10                 |
| In-Service Cover                            | 76901-40                 |
| Draft Shield Snap Clamp (TS120S and TS400D) | 77334-00                 |
|   |                          |

#### **ACCESSORIES**

| Draft Shield Kit<br>Chamber Size: 6.0"W x 6.375"L x 4.25"H | OHAUS<br><u>Part No.</u><br>76934-01                                 |
|--|--|
| RS-232 Interface Kit                                       | 77018-01   |
| Cable with PRINT switch for RS232 Interface                | AS017-26   |
| Leveling Foot Conversion Kit:                              | 77327-00   |
| Weigh Below Hook   | 76790-00   |
| In-Service Cover Kit                                       | 76901-30   |
| Anti Theft Device  | 76288-00   |
| Calibration Weights: 40g 50g 100g 200g 400g 2kg 4kg        | 49044-01<br>49054-01<br>49015-01<br>49025-01<br>49045-01<br>49026-01 |

#### LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.



## PRECISION Standard Electronic Balances TS Series

### **Instruction Manual**

**NOTE:** THIS EQUIPMENT HAS BEEN TESTED AND FOUND TO COMPLY WITH THE LIMITS FOR A CLASS A DIGITAL DEVICE, PURSUANT TO PART 15 OF THE FCC RULES.

THESE LIMITS ARE DESIGNED TO PROVIDE REASONABLE PROTECTION AGAINST HARMFUL INTERFERENCE WHEN THE EQUIPMENT IS OPERATED IN A COMMERCIAL ENVIRONMENT. THIS EQUIPMENT GENERATES, USES, AND CAN RADIATE RADIO FREQUENCY ENERGY AND, IF NOT INSTALLED AND USED IN ACCORDANCE WITH THE INSTRUCTION MANUAL, MAY CAUSE HARMFUL INTERFERENCE TO RADIO COMMUNICATIONS. OPERATION OF THIS EQUIPMENT IN A RESIDENTIAL AREA IS LIKELY TO CAUSE HARMFUL INTERFERENCE IN WHICH CASE THE USER WILL BE REQUIRED TO CORRECT THE INTERFERENCE AT HIS OWN EXPENSE.

THIS DIGITAL APPARATUS DOES NOT EXCEED THE CLASS A LIMITS FOR RADIO NOISE EMISSIONS FROM DIGITAL APPARATUS AS SET OUT IN THE INTERFERENCE-CAUSING EQUIPMENT STANDARD ENTITLED "DIGITAL APPARATUS", ICES-003 OF THE DEPARTMENT OF COMMUNICATIONS.

CET APPAREIL NUMERIQUE RESPECTE LES LIMITES DE BRUITS RADIOELECTRIQUES APPLICABLES AUX APPAREILS NUMERIQUES DE CLASSE A PRESCRITES DANS LA NORME SUR LE MATERIEL BROUILLEUR: "APPAREILS NUMERIQUES", NMB-003 EDICTEE PAR LE MINISTRE DES COMMUNICATIONS.

Unauthorized changes or modifications to this equipment are not permitted.

#### **TABLE OF CONTENTS**

| INT | RODUCTION                           | 5  |
|-----|-------------------------------------|----|
| DES | SCRIPTION                           | 5  |
| UNI | PACKING                             | 6  |
| INS | TALLATION                           | 7  |
|     | Environment                         | 7  |
|     | Draft Shield                        | 7  |
|     | Pan and Pan Support                 | 8  |
|     | AC Adapter                          | 8  |
| OPI | ERATION                             | 9  |
|     | Turning the Balance ON              | 9  |
|     | Stabilization                       | 9  |
|     | Auto Range Models                   | 9  |
|     | Checking Calibration                | 10 |
|     | Weighing 1                          | 10 |
|     | Taring 1                            | 11 |
| USI | NG MENUS TO CONFIGURE THE BALANCE 1 | 12 |
| CAI | LIBRATION MENU 1                    | 13 |
|     | Calibration Menu Protection         | 13 |
|     | Calibration Weights                 | 13 |
|     | Span Calibration                    | 14 |
|     | Linearity Calibration 1             | 15 |
|     | End 1                               | 16 |
| USI | ER MENU 1                           | 17 |
|     | User Menu Protection                | 17 |
|     | Reset to Factory Defaults           | 18 |
|     | Averaging Level 1                   | 19 |
|     | Stability Range                     | 20 |
|     | Auto-Zero                           | 20 |
|     | End                                 | 21 |


| SETUP MENU               | 22 |
|--------------------------|----|
| Setup Menu Protection    | 22 |
| Unit Selection           | 23 |
| Lock Switch              | 24 |
| End                      | 25 |
| MENU LOCK-OUT PROTECTION | 25 |
| CARE AND MAINTENANCE     | 26 |
| TROUBLESHOOTING          | 26 |
| Error Codes              | 27 |
| SERVICE INFORMATION      | 28 |
| SPECIFICATIONS           | 29 |
| PARTS INFORMATION        | 30 |
| REPLACEMENT PARTS        | 30 |
| ACCESSORIES              | 30 |
| LIMITED WARRANTY         | 31 |

#### INTRODUCTION

This manual covers installation, operation and troubleshooting for the Ohaus Precision Standard balances, Models TS120, TS200, TS400, TS400D, TS600, TS2K, TS4K and TS4KD. To insure proper operation of the balance, please read this manual completely.

#### DESCRIPTION

The Ohaus Precision Standard series balances are precision weighing instruments, designed to provide years of service with virtually no maintenance. The Precision Standard series is constructed using a die-cast aluminum base finished with a durable epoxy powder paint which is resistant to commonly used acids, contains a one piece solid-state precision electronics PC board, a seven digit LCD display which is 0.6 inches in height. All Precision Standard series balances are factory set to measure in grams. Each balance operates through a series of menus which enable precise calibration and linearity along with various other parameters which enhances operation. A built in lock switch prevents preset settings from being changed. To prevent measurements from being affected by air currents, a draft shield is used on Models TS120, TS200 and TS400D balances. Power is supplied through an AC adapter which is available in five voltages for world-wide usage. Accessories include: an RS232 interface kit which allows printing of results through an external computer, an RS232 Interface cable with a print switch, weigh below hook, security device and calibration weights.

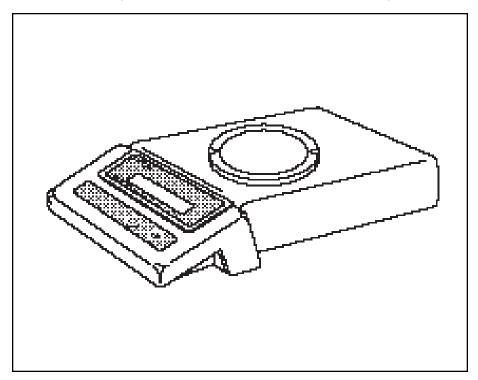


PRECISION Standard Balance

#### **UNPACKING**

Your Precision Standard balance was shipped with the following items:

- a pan
- · a pan support
- an AC power adapter
- a draft shield (TS120, TS200 and TS400D only) includes draft shield and snap clamp
- · this instruction manual
- · your warranty card

It is recommended to save the carton and packing material for storing, transporting the balance or returning it for service.

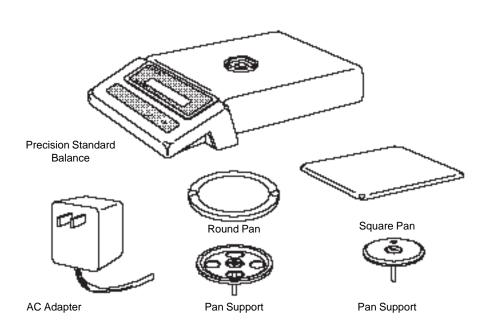

(Round pan and support for Models TS120, TS200, TS400, TS600, and TS400D) (Square pan and support for Models TS2K, TS4K and TS4KD)

#### **INSTALLATION**

#### **Environment**

The balance should always be used in an environment which is free from excessive air currents, corrosives, vibration, and temperature or humidity extremes. These factors will affect displayed weight readings.

#### DO NOT install the balance:

- Next to open windows or doors causing drafts or rapid temperature changes.
- Near air conditioning or heat vents.
- · Near vibrating, rotating or reciprocating equipment.
- · Near magnetic fields or equipment that generates magnetic fields.
- On an unlevel work surface.


#### Draft Shield (TS120, TS200 and TS400D)

To install the draft shield:

 Position the draft shield on top of the balance as shown. Make sure the rib at the front of the draft shield base butts up against the raised lip of the display panel.

2. The snap clamp should be oriented so that the double clip is toward the front of the balance.

Insert the double clip of the snap clamp into the opening in the draft shield base, and clip the draft shield base to the balance. Press the rear clip of the snap clamp into place.

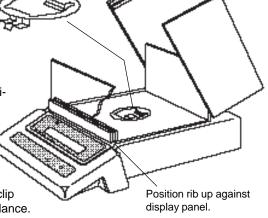


#### **Pan and Pan Support**

#### Square Pan

Insert the pan support into the hole in the weighing mechanism as shown in the illustration. Make sure the hole in the pan support faces the rear of the balance. Once installed, the pan support should not rotate.


The pan has a guide pin which protrudes from the bottom. Place the pan on the support making sure the guide pin is inserted in the hole in the pan support.

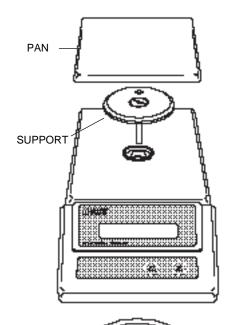

#### **Round Pan**


Insert the pan support into the hole in the weighing mechanism.

Place the pan on the support.

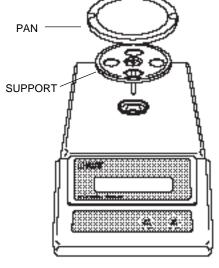

#### **AC Adapter**

Plug the molded connector of the adapter into the receptacle at the rear of the balance. Plug the adapter into a convenient AC outlet. When power is applied to the balance, it will begin a self test cycle. During this time, the display will count downfrom 10 and display the word CHEC.

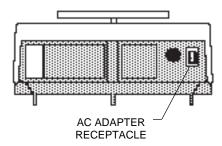


#### **OPERATION**

#### **Turning the Balance ON**

With no load on the pan, turn the balance ON by pressing the ON TARE button. When first switched ON, all segments of the display should be on as shown in the illustration.

This display check will be displayed briefly, then the model number of the balance, followed by a short countdown. The display will momentarily blank and then indicate zero. All of the displays shown at the right only appear during initial power-up.



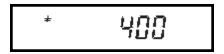




•

#### **Stabilization**

Before initally using the balance, allow time for it to adjust to changes in environment. The balance need only be plugged in to warm up. Recommended warm up period is thirty minutes.

## C 33K3

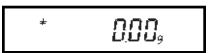

## Auto Range Models (TS400D and TS4KD)

Auto range balances offer both a fine range (lower capacity/higher readability) and a coarse range (higher capacity/lower readability). When first turned on, the balance is in the fine range. It remains in this range until the weight on the pan exceeds the fine range capacity. When weight on the pan is greater than the fine range capacity, the balance switches to the coarse range.

If weight on the pan falls back to within the fine range capacity, coarse range readability remains in effect until you tare the balance with no weight on the pan.


#### **Checking Calibration**

Before using the balance, calibration should be checked. The balance has been calibrated before shipment, however, it could be influenced by factors such as:

- Variations in the earths gravitational field at different latitudes of the world.
- · Rough handling.
- Changes in work location.
- · Height above sea level.


To check the balance's calibration, place a known mass on the center of the pan and read the displayed weight.

If the displayed weight differs from the known weight by more than acceptable limits, refer to the Calibration Menu and the Specifications at the rear of the manual.

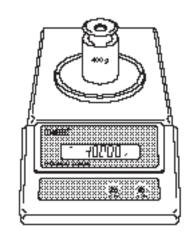

#### Weighing

- 1. Press ON to rezero the display.
- 2. Place the object(s) or material to be weighed on the pan.
- 3. Wait for the stability indicator to appear before reading the weight.

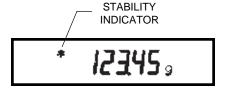

#### **Taring**

When weighing material or objects that must be held in a container, taring enables you to store the container weight in the balance's memory, separate from the weight of the material in the container.


- Place an empty container on the pan. Its weight will be displayed.
- 2. Press ON TARE


The display will show zero and the container's weight will be stored in memory.

- Add material to the container. As material is added, its net weight will be displayed.
- Removing the container and material from the pan will cause the balance to display the container's weight as a negative number.

Tared weight remains in balance memory

until (ON tare) is pressed again.

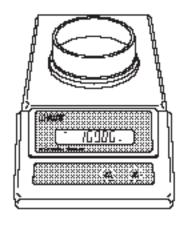

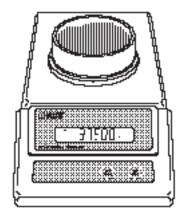

#### **USING MENUS TO CONFIGURE THE BALANCE**


Precision Standard balances contain four display menus which enable you to calibrate and configure the balance for your specific operating requirements.

**Calibration Menu:** Used to calibrate the balance for span or linearity.

**User Menu**: Used to adapt balance to environmental conditions.

**Setup Menu:** Used to enable, disable or customize different balance features.

**Print Menu:** Used to configure the RS232 interface.

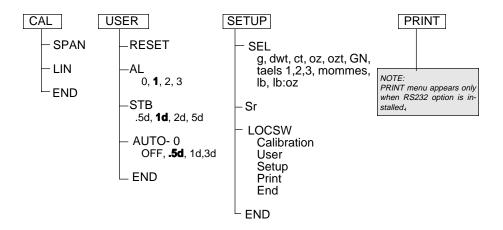

To access a menu, press and hold  $\binom{\text{ON}}{\text{TARE}}$  until desired menu appears, then release it.

Original factory default settings are shown in boldface type.

Use these buttons to step through menus and select submenus:



#### **CALIBRATION MENU**

Precision Standard balances can be calibrated in two ways: Span calibration or Linearity calibration. Span calibration resets the balance's weighing range using two weight values: zero and a weight value at or near the balance's capacity. Linearity calibration minimizes deviation between actual and displayed weights within the balance's weighing range. Three weight values are used: zero, a weight value within the balances weighing range, and a weight value at or near the balance's specified capacity. The following table shows the sequence in which submenus appear on the Calibration menu.

| CALIBRA | ATION | MENU | <b>TABLE</b> |
|---------|-------|------|--------------|
|---------|-------|------|--------------|

| SPAn | Selects span calibration.          |
|------|------------------------------------|
| Lin  | Selects linearity calibration.     |
| End  | Used to exit the Calibration menu. |

#### **Calibration Menu Protection**

The calibration menu may be locked out to prevent unauthorized personnel from changing calibration. To lock out the calibration menu, refer to section titled Menu Lock-Out Protection.

**NOTE**: If calibration has been locked out, you will not be able to access it.

#### **Calibration Weights**


Before beginning calibration, make sure weights are on hand. If you begin calibration and realize weights are not available, either turn the balance off, or go through the procedure without weights. The balance will use previously stored calibration data. Calibration should be performed as necessary to ensure accurate weighing. Weights required to perform the procedures are listed in the adjacent table.

| <b>CALIBRATION WEIGHTS</b> |                      |                     |  |
|----------------------------|----------------------|---------------------|--|
| MODEL                      | LINEARITY<br>WEIGHTS | SPAN ONLY<br>WEIGHT |  |
| TS120                      | 50g, 100g            | 100g                |  |
| TS200                      | 100g, 200g           | 200g                |  |
| TS400                      | 200g, 400g           | 400g                |  |
| TS400D                     | 50g, 400g            | 400g                |  |
| TS600                      | 200g, 500g           | 500g                |  |
| TS4K                       | 2kg, 4kg             | 4kg                 |  |
| TS2K                       | 1kg, 2kg             | 2kg                 |  |
| TS4KD                      | 500g, 4kg            | 4kg                 |  |

It is recommended that masses must meet or exceed ASTM Class 1 Tolerance. Calibration masses are available as accessories.

#### **Span Calibration**


Press and hold (TARE) until CAL is displayed, then release it. Balance will display SPAN.

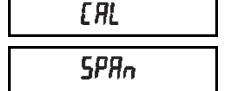



- 2. Press ON TARE.
- When (TARE) is released, C 0g will be displayed indicating that no weight should be on the pan.
- 4. Press (TARE). The display will show
   -C-followed by the value of the weight which must be placed on the pan.

NOTE: Do not disturb the balance when -C- is displayed. Disturbances will result in improper calibration.

- Place the required weight on the pan and press ON TARE. The display will show -C- while the balance recalibrates.
- When the weight on the pan is displayed along with the current unit indicator, the balance is recalibrated.

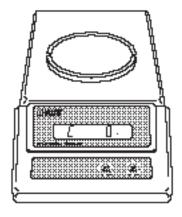

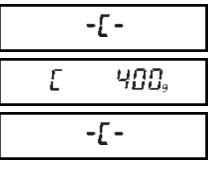


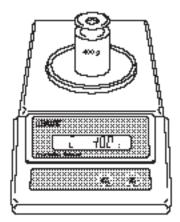


#### **Linearity Calibration**


- Turn Balance on. After zero reading, press and hold ON UNTARE until CAL is displayed, then release. Balance will display SPAN.
- 2. Press OFF and the display will show LIN.
- 3. Press TARE to start the Linearity Calibration Procedure. When the is released, C 0g will be displayed, indicating that no weight should be in the pan.
- 4. Press (NAME). The display will show
   -C-followed by the value of the weight which must be placed on the pan.
- 5. Place the required weight on the pan.

EAL SPAn L in

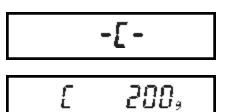

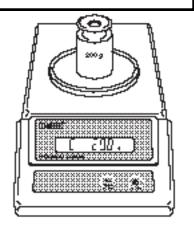



6. Press ON TARE. The display will show -C- momentarily, then C followed by the next weight to be placed on the pan. Do not disturb the balance when -C- is displayed. Disturbances will result in improper calibration.

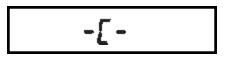

E 400,

7. Place the required weight on the pan, then press ON TARE. The display will show -C- while the balance recalibrates. When the weight on the pan is displayed along with the current indicator, the balance is recalibrated.

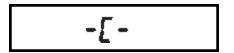

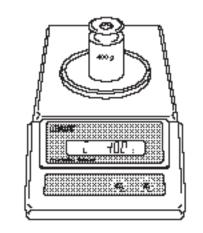

#### **End**

If you have entered the Calibration menu and do not wish to calibrate the balance, use END to return to normal weighing operations.

Repeately press wode until End is displayed.

Press (ON ), when released, the balance

will returned to normal weighing opera-

End

#### **USER MENU**

The User menu is used to adapt the balance to environmental conditions. It contains submenus which enable you to reset the balance to factory default settings or to select specific range settings. Access to the User menu can be disabled using the Lock Out switch. The following table shows the sequence in which submenus appear on the User menu.

#### **USER MENU TABLE**


| rESEŁ  | Sets all submenus below to original factory default settings. Reset does not appear if menu has been locked out. |
|--------|--|
| AL     | Specifies the averaging level.   |
| 269    | Specifies the desired stability range.   |
| Auto-0 | Sets Auto-Zero threshold.  |
| End    | Used to exit the Setup menu and store your selections.   |

#### **User Menu Protection**


The User menu may be locked out to prevent unauthorized personnel from changing the settings. To lock out the User menu, refer to the section titled Menu Lock-Out Protection.

USEr

**NOTE**: If -SAFE- is displayed, the User menu has been locked out. Settings may be viewed but not changed. See the Menu Lock-Out Protection section to enable it for making changes.

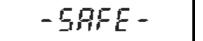

To access the User menu, press and hold

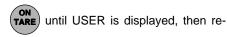

lease it.

To access a submenu:

Repeatedly press (NOTE) until the desired submenu is displayed.

Press ON to select the displayed submenu.

**NOTE**: You must use END to store any changes you make to the User menu.

The following sections describe each item on the User menu in detail.

#### **Reset to Factory Defaults**

This submenu enables you to reset all User menu selections to the factory default settings outlined in the adjacent table.

To reset to factory defaults:


- 1. Access the RESET submenu.
- 2. Press MODE to change the setting.

Select YES to reset settings or, no to leave current settings.

Press ON to accept the displayed setting.

rESEŁ

| USER MENU<br>FACTORY DEFAULTS |      |  |
|-------------------------------|------|--|
| Averaging Level               | AL 1 |  |
| Stability Range               | 1d   |  |
| Auto-Zero Tracking            | .5d  |  |

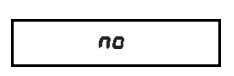

## **Averaging Level**

Averaging level compensates for vibration or excessive air currents. During operation, the balance continually takes weight readings from the weighing cell. Successive readings are then digitally processed to achieve a stabilized display. Use this submenu to specify how much processing you need to obtain stable results.

NOTE: Averaging level does not affect balance accuracy.

Select one of four averaging levels using the adjacent table as a guide.

To view or change the averaging level:

- 1. Access the AL submenu to display the current setting.
- 2. Press MODE to change the setting.

Press ON to accept the displayed setting.

When on is released, AL will be displayed again and the Setup menu will be returned.

## AL

#### **AVERAGING LEVEL**

AL 0 reduced stability, fastest stabilization time

## AL 1 normal stability, normal stabilization time

- AL 2 more stability, slow stabilization time
- AL 3 maximum stability, slowest stabilization time



## **Stability Range**

The stability range specifies how much a displayed weight may change while the stability indicator remains ON. When displayed weight changes beyond the allowable range, the stability indicator turns OFF indicating an unstable condition. Precision Standard balances permit you to select one of four stability ranges (in divisions) as shown in the table.

When the RS232 interface is configured to print stable data only, the stability range also governs data output. Displayed data will only be output if it is within the selected stability range.

To view or change the stability range:

- 1. Access the Stb submenu to display the current setting.
- 2. Press OFF to change the setting.

Press ON to accept the displayed setting.

When TARE is released, Stb will be displayed again and the Setup menu will be returned.

#### **Auto-Zero**

Auto-Zero minimizes the effects of temperature changes and shift on the zero reading. By defining a threshold level in divisions, the balance maintains the zero display until the threshold is exceeded. This submenu permits you to select one of three threshold levels, or turn the feature OFF. Auto-Zero only functions when the display reads zero.

SEB

#### **STABILITY RANGE**

.5d smallest range: stability indicator is ON only when displayed weight is within .5 divisions

## 1d reduced range

- 2d normal range
- 5d largest range: stability indicator is ON even though displayed weight changes slightly

1 8

Auto-0

To view or change the Auto-Zero setting:

 Access the Auto-0 submenu to display the current setting.

5 d

2. Press OFF to change the setting.

Press ON to accept the displayed setting.

When TARE is released, Auto-0 will be displayed again and the User menu will be returned.

#### **AUTO ZERO**

OFF turns Auto-Zero OFF

.5d sets threshold to .5 divisions1d sets threshold to 1 division

3d sets threshold to 3 divisions

### End

You must use END to exit the User menu. Changes you make in the User menu are only stored in memory if you use END.

To exit the User menu and store your settings, press (ON ) when End is displayed.

When ON is released, the balance will be returned to normal weighing operations.

End

#### **SETUP MENU**

The Setup menu enables you to retain program balance parameters once they have been set. Access to the Setup menu can be disabled using the Lock Out switch. The following table shows the sequence in which submenus appear on the Setup menu.

#### **SETUP MENU TABLE**

| CCI         | Enables the selection of weighing     |
|-------------|---------------------------------------|
| 200         | units.                                |
| 5           | Indicates the software version of the |
|             | balance.                              |
| L 🛮 🖰 ԵՐԱ   | Enables individual or all menus to be |
| F D F 2 D D | locked out.                           |
| C _ J       | Used to exit the Setup menu and       |
| End         | store your selections.                |

## **Setup Menu Protection**

The Setup menu may be locked-out to prevent unauthorized personnel from changing settings. To lock out the Setup menu, refer to the section titled Menu Lock-Out Protection.

**NOTE**: If -SAFE- is displayed, the Setup menu has been locked out. Settings may be viewed but not changed. See the Menu Lock-Out Protection section to enable it for making changes.

To access the Setup menu, press and hold ON until SETUP is displayed, then

**-**58EUP-

release it.

To access a submenu:

-5888-

- Repeatedly press (MODE) until the desired submenu is displayed.
- 2. Press (ON to select the displayed

submenu.

**NOTE**: You must use END to store any changes you make to the Setup menu.

The following sections describe each item on the Setup menu in detail.

#### **Unit Selection**


Unit selection permits you to specify which weighing units will be enabled for use during operation. The adjacent table lists the units available.

To enable or disable the various weighing units, use the following procedure:

- 1. Access the SEL menu.
- The display will show the grams unit indicator (g) along with the current status (ON/OFF).
- 3. Press MODE to change the status.
- 4. Press (ON TARE) to accept the displayed status. When (ON TARE) is released, the display will show the next unit indicator with the current status.
- 5. Set each unit ON or OFF as in step 3.

# SEL

| Weighing Units   |           |  |
|------------------|-----------|--|
| g grams : mommes |           |  |
| dwt pennyweight  | ct carats |  |
| oz t troy ounces | t taels   |  |
| lb               | lb:oz     |  |
| gn               | OZ        |  |

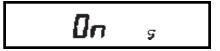



#### **Taels**

If taels are enabled, you will be required to choose one of three different taels: Hong Kong, Singapore, or Taiwan.

When the display shows TAEL 1, press




to change to another tael, press



to accept the displayed tael.

When the last weighing unit has been set, the display will show SEL again and the Setup menu will be returned.

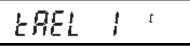

(Hong Kong)

FWET 5 .

(Singapore)

F867 3 .

(Taiwan)

#### Lockswitch

Lockswitch enables you to lock out one or more menu selections. Each menu can be individually locked on or off after all functions have been set. The Calibration, User, and Setup menus can be individually locked on or off by selecting the appropriate menu and then locked by the switch located under the right hand side of the control panel. Before performing the lock out procedure, decide which functions of the balance are to be locked on or off.

LOCSUJ

- Access the LOCSW submenu. When ON TARE is released, the LOCSW submenu is displayed.
- To access one or more menus, press
   ON TARE to select the calibration menu,
   -CAL- is displayed.

**NOTE**: Pressing MODE changes the selection to the other menus.

3. To select a YES or NO, press MODE

**NOTE**: The MODE switch acts as a toggle and can select either YES or NO.

To confirm your selection, press TARE
 again. The display indicates the last menu you were in.

455

no

5. To lock out the other menus, press



and repeat the procedure in

steps 3 and 4.

#### **End**

You must use END to exit the Setup menu. Changes you make in the Setup menu are only stored in memory if you use END.

To exit the Setup menu and store your settings, press ON when END is displayed.


When tiare is released, the balance will be returned to normal weighing operations.

## MENU LOCK-OUT PROTECTION

Access to the Calibration, User, and Setup menus, can be disabled using the lock out switch located under the right side of the balance, near the display.

- 1. Turn the display off and unplug the power cord.
- Slide the balance toward you, with the front over the edge of a table. (You can also turn the balance on its left side, but if you do, you MUST remove the pan and spill ring first!)
- Locate hole under display where switch is located.
- Using a small screwdriver, slide the switch forward for LOCKED or back for UNLOCKED.
- 5. Plug in the power cord and turn on the balance.

## End

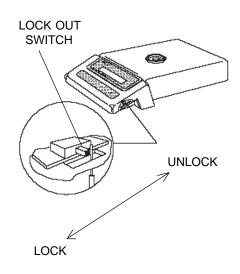

## **CARE AND MAINTENANCE**

To keep the balance operating properly, the housing and pan should be kept clean and free from foreign material. If necessary, a cloth dampened with a mild detergent may be used. Keep calibration weights in a safe dry place.

## **TROUBLESHOOTING**

| SYMPTOM  | PROBABLE CAUSE(S)  | REMEDY  |
|--|--|---|
| Unit will not turn on.                         | Power adapter not plugged in or properly connected to balance. | Check power adapter connections.  |
| Incorrect weight reading.                      | Balance was not re-<br>zeroed before weighing.                 | Press ON TARE with no weight on the pan, then weigh item.                               |
|  | Balance not properly calibrated.                               | Recalibrate correctly.  |
| Unable to store menu settings/changes.         | END not being used to exit menus.                              | You MUST use END to exit menus and save settings.                                       |
| RS232 interface not working (when installed)   | Print menu settings not properly set up.                       | Verify interface settings in Print menu correspond to those of peripheral device.       |
|  | Cable connections.   | Check cable connections.  |
| Random segments displayed or display locks up. | Microprocessor lock up.  | Unplug the power cord, then replug again. If condition persists, unit must be serviced. |
| Unable to change settings.                     | Lock set ON.   | Set Lock switch to OFF.   |
| Unstable readings.                             | Vibration on table surface.                                    | Place balance on a stable surface or change averaging level.                            |
| Error message display.                         |  | See Error Codes Table.  |

#### **Error Codes**

The following list describes the various error codes and which can appear on the display and the suggested remedy.

#### **Data Errors**

0.0 Transient error (hardware error, probably static discharge). If error persists, the balance must be serviced.

#### **Tare Errors**

2.0 Balance is unable to stabilize within time limit after taring. Environment is too hostile or balance needs recalibration.

#### **Calibration Errors**

3.0 Incorrect or no calibration weight used for calibration. Recalibrate with correct weights.

#### **RS232 Errors**

- 4.0 Bad RS232 frame. Check RS232 menu parameters and correct.
- 4.4 RS232 buffer is full (if installed). May occur if no printer or computer is connected to the interface. To clear buffer, turn balance off or enter Print menu and select END.
- 4.5 Function is disabled by the Lock switch.

#### **User Errors**

7.2 Number outside of display capacity.

#### **Over-Under Load Errors**

- 8.0 Hardware error causing an internal weight signal which is too low. Check if pan or pan support is off. If not, the balance must be serviced.
- 8.1 Hardware error caused by an internal weight signal which is too high. Check load on the platform which may be excessive. If error persists, the balance must be serviced.
- 8.2 Power-on load out of specification: Balance was turned on with load on pan or pan off balance. No load may be on pan when turned on and pan must be in place.
- 8.3 Rated capacity exceeded. Remove excessive weight from pan.
- 8.4 Underload condition on balance. Check that the proper pan and pan support are installed.

## **Error Codes (Cont.)**

#### **CheckSum Errors**

- 9.0 Bad factory checksum. If error persists, have the balance serviced.
- 9.5 Bad factory calibration checksum. If error persists, have the balance serviced.
- 9.6 Bad mode checksum. Turn the balance off using the front panel controls. If the error persists, have the balance serviced.
- 9.7 Invalid setup data checksum. Check Setup, User and Print menu (when RS232 is installed) settings. If possible, try to enter menus and exit using END to restore menu settings. May be caused by a faulty component, or in rare cases, a severe static charge. If error persists, balance must be serviced.
- 9.8 Hardware error causing invalid calibration data checksum. Balance may need recalibration - particularly linearity calibration. If error persists, balance must be serviced.
- 9.9 Invalid temperature compensation checksum. Balance will work with default temperature compensation data, however, error will occur each time balance is turned on. Have balance serviced.

## **SERVICE INFORMATION**

If the Troubleshooting section does not resolve or describe your problem, you will need to contact an authorized Ohaus Service Agent. For Service assistance in the United States, please call Ohaus Corporation toll-free at (800) 526-0659, an Ohaus Product Service Specialist will be available to help you.

| MODEL                             | TS120   | TS200  | TS400D               | TS400      | TS600        | TS2K     | TS4KD     | TS4K |
|-----------------------------------|---|--------|----------------------|------------|--------------|----------|-----------|------|
| Capacity (g)                      | 120   | 200    | 400/80               | 400        | 600          | 2000     | 4000/800  | 4000 |
| Readability (g)                   | 0.0   | 01     | 0.01/0.001           |            | 0.01         |          | 0.1/0.01  | 0.1  |
| Weighing mode                     |   | g, Ib  | o, oz, lb:oz, c      | t, dwt, ta | els, oz t, ç | gn, momn | ne        |      |
| Repeatability<br>(Std. dev.) (g)  | 0.00  | )1     | 0.007/0.001          | 0.0        | 007          | 0.01     | 0.07/0.01 | 0.07 |
| Linearity (g)                     | ±0.001  | ±0.002 | ±0.01/0.001          | ±0         | ).01         | ±0.02    | ±0.1/0.01 | ±0.1 |
| Tare range                        |   |        | Full ca <sub>l</sub> | oacity by  | subtraction  | on       |           |      |
| Stabilization time                |   |        | 2.!                  | second     | S            |          |           |      |
| Sensitivity drift (10 - 30 °C)    | 10ppm/ °C   |        |                      |            |              |          |           |      |
| Operating temperature             | 50° to 104°F/10 ° to 40°C                               |        |                      |            |              |          |           |      |
| Calibration                       | External digital calibration                            |        |                      |            |              |          |           |      |
| Display (in/cm)                   | LCD (0.6/1.5 high)                                      |        |                      |            |              |          |           |      |
| Power requirements                | AC adapter: 100, 120, 220, 240 V ac, 50/60 Hz           |        |                      |            |              |          |           |      |
| Platform size (WxHxD) (in/cm)     | 4.8/ 6 x 6.5/<br>12.2 15.2 x 16.5                       |        |                      |            | 5            |          |           |      |
| Dimensions<br>(W x H x D) (in/cm) | 8 x 3.75 x 14/20.3 x 9.5 x 35.6<br>without draft shield |        |                      |            |              |          |           |      |
| Net weight (lb/kg)                | 9.8/4.4   |        |                      |            |              |          |           |      |
| Shipping weight (lb/kg)           | 13.6/6.2  |        |                      |            |              |          |           |      |

## **PARTS INFORMATION**

If you require replacement parts or would like to purchase accessories, please call Ohaus Corporation toll-free at (800) 526-0659, an Ohaus Product Parts Specialist will be available to help you.

## **REPLACEMENT PARTS**

| Description AC Adapters:                          | OHAUS<br><u>Part No.</u> |
|---|--------------------------|
| 100V  | 90766-01                 |
| 120V  | 90765-01                 |
| 220V  | 90767-01                 |
| 240V  | 90768-01                 |
| 240V Australia                                    | 90524-15                 |
| Pan - 4.7" dia.                                   | 77262-10                 |
| Pan - 6.0" x 6.5"                                 | 77298-10                 |
| Draft Shield Snap Clamp (TS120, TS200 and TS400D) | 77334-00                 |

## **ACCESSORIES**

| In-Service Cover   | OHAUS<br><u>Part No.</u><br>78211-02                                 |
|--|--|
| Draft Shield Kit<br>Chamber Size: 6.0"W x 4.25"H x 6.375"D                   | 76934-03   |
| RS232 Interface Kit  | 77018-01   |
| Cable for RS232 Interface  | AS020-17   |
| Cable with PRINT switch for RS232 Interface                                  | AS017-25   |
| Leveling Foot Conversion Kit:  | 77327-00   |
| Weigh Below Hook   | 76790-00   |
| Security Device  | 76288-01   |
| Calibration Masses - ASTM Class 1 Tolerance: 50g 100g 200g 400g 500g 1kg 2kg | 49054-11<br>49015-11<br>49025-11<br>49045-11<br>49055-11<br>49016-11 |
| 4kg  | 49046-11   |

#### LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.



Ohaus Corporation 29 Hanover Road, Florham Park, NJ 07932, USA

Tel: (973) 377-9000, Fax: (973) 593-0359

With offices worldwide.



# Weigh Below Hook Accessory


P/N 76790-00

For PRECISION *Standard*Flectronic Balances

**Installation Instructions** 

## INSTALLATION

- 1. Turn the balance OFF.
- Remove the pan.
- 3. Hold the pan support in place and turn the balance over or on its side.
- 4. Screw the hook into the threaded hole in the pan support which is visible through the access hole in the bottom of the balance.

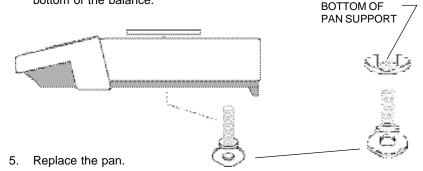

OHAUS® is the registered trademark of Ohaus Corporation as are the following trademarks: AUTOGRAM®, BRAINWEIGH®, CENT-O-GRAM®, CHECK-O-GRAM®, CUBE-O-GRAM®, DEC-O-GRAM®, DIAL-O-GRAIN®, DIAL-O-GRAM®, DU-O-MEASURE®, GALAXY®, LUME-O-GRAM®, PORT-O-COUNT®, PORT-O-GRAM®, PRACT-O-GRAM®, PRIMER®, STO-A-WEIGH®, TOUCH-N-WEIGH®, 5-0-5®, and 10-10®.



#### **OHAUS CORPORATION**

P.O. Box 900 29 Hanover Road Florham Park, N.J. 07932 Tel: 201-377-9000 Telex: 6853191 OHAUS UW

Fax: 201-593-0359

#### **OHAUS EUROPE, LTD.**

Broad Lane, Cottenham Cambridge CB4 4SW ENGLAND Tel: 0954-51343 Telex: 817285 OSCALE G Fax: 0954-50205



# Draft Shield Accessory Kit

P/N 76934-01

For PRECISION *Standard* and *Plus*Electronic Balances

Models TS400, TP400 and TP600

**Installation Instructions** 

### **UNPACKING**

Verify that a Draft Shield and Snap Clamp have been included in the kit.

INSTALLATION Double Clip 1. Remove the pan and pan support. Snap Position the draft shield Clamp on top of the balance as shown. Make sure the rib at the front of the draft shield base butts up against the raised lip of the display panel. 3. The snap clamp should be oriented so that the double clip is toward the front of the balance. Cover First, insert the double clip Position rib up against end of the snap clamp fully into display panel. the opening towards the front of the draft shield base and the underside of the cover. Next, press the rear clip of the

4. Replace the pan and pan support.

snap clamp down into place.

OHAUS $^{\circ}$  is the registered trademark of Ohaus Corporation as are the following trademarks: CENT-O-GRAM $^{\circ}$ , CHECK-O-GRAM $^{\circ}$ , DEC-O-GRAM $^{\circ}$ , DIAL-O-GRAM $^{\circ}$ , PRIMER $^{\circ}$ , 5-0-5 $^{\circ}$ , and 10-10 $^{\circ}$ .

#### **OHAUS CORPORATION**

P.O. Box 900 29 Hanover Road Florham Park, N.J. 07932, USA Tel: 201-377-9000 Telex: 6853191 OHAUS UW Fax: 201-593-0359

With other offices in England, Mexico, Canada, Japan, Germany, France, Switzerland





# Leveling Kit Accessory

P/N 77327-00

# For PRECISION Standard Electronic Balances

## Installation Instructions

## **UNPACKING**

The Leveling Kit includes the following components:

1 FCC label 2 lockwashers 1 level assembly with 3 set screws installed

2 phillips screws


1 brass foot

4 leveling feet (2 pairs)

## INSTALLATION


A phillips screwdriver, 1/16" hex wrench and 9/64" hex wrench will be required to install the leveling kit.

- 1. Disconnect power to the balance.
- 2. Remove the pan and pan support.
- Turn the balance over.
- Remove the FCC label. The label covers two clearance holes that will be used for mounting the leveling assembly.
- 5. Peel off the two rubber feet at the front of the balance.
- Using a 9/64" hex wrench, remove the socket head screw at the rear of the balance. Leave the washer in place.

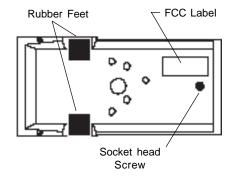


- Install the brass foot in place of the socket head screw.
- Screw the leveling feet in leaving approximately 3/16" clearance between the bottom of the balance and the foot.

NOTE: Two pairs of leveling feet are included in this kit. The all plastic feet are for earlier TS Models. The leveling feet with metal threaded shafts are for later versions.

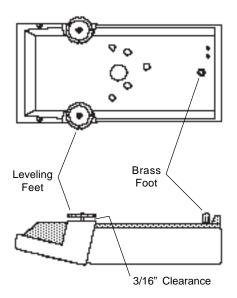


- Remove the three phillips head screws which secure the balance cover to the base.
- 10. Hold the base and cover together and turn the balance right side up.
- 11. Lift the cover off.

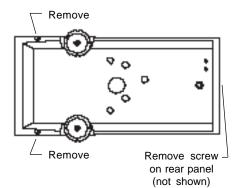

Lockwashers-

12. Using the two phillips screws and lockwashers provided in the kit, fasten the leveling assembly as shown in the diagram.


Screws should be snug but not tight. DO NOT COMPLETELY TIGHTEN THE SCREWS.

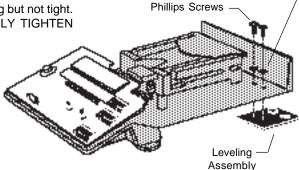

- 13. Place the balance on a level work surface. ENSURE SURFACE IS LEVEL BEFORE CONTINUING. Position the balance so that the leveling assembly extends from the edge of the work surface.
- 14. Using a 1/16" hex wrench adjust the three set screws in the leveling assembly as needed to center the bubble in the level indicator.

When the bubble is centered, tighten the two phillips screws.

 Apply the new FCC label (included in the kit) to the bottom of the balance.
 Do not to cover any screws or holes.

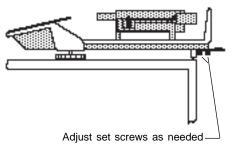



- Replace the cover and fasten with the three phillips screws that were removed in step 9.
- 17. Replace the pan support and pan.
- 18. Place the balance in the location where it will be used and verify it is level. If minor adjustment is needed, use the leveling feet.
- Recalibrate the balance as explained in the instruction manual supplied with the balance.

Ohaus Corporation 29 Hanover Road, Florham Park, NJ 07932, USA

Tel: (973) 377-9000, Fax: (973) 593-0359

With offices in England, Germany, France, Spain, Italy, Poland, Mexico, Japan, Korea and China.



# In-Service Cover Accessory Kit

P/N 76901-30

For PRECISION *Standard* and *Plus* Electronic Balances

## **Installation Instructions**




## **UNPACKING**

The kit includes a cover and one adhesive pad.

## **INSTALLATION**

- 1. Remove the pan and pan support from the balance.
- Place the cover over the balance as shown in the diagram. If desired, the adhesive pad may be used to secure the cover to the balance to avoid possible movement.

Peel the backing off of one side of the adhesive pad, center adhesive pad over hole in top cover and press the pad onto the top of the balance as shown in the diagram.

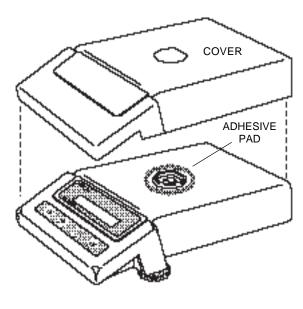

Remove the backing from the top of the pad, place the cover on the balance and press down on the pad.

3. Replace the pan and pan support.

Ohaus Corporation 29 Hanover Road, Florham Park, NJ 07932, USA

Tel: (973) 377-9000, Fax: (973) 593-0359

With offices in England, Germany, France, Spain, Italy, Poland, Mexico, Japan, Korea and China.



RS232 INTERFACE ACCESSORY
Part Number 77018-01
For
ANALYTICAL Standard Series
and
PRECISION Standard Balance Series

## **Instruction Manual**

**NOTICE:** THIS EQUIPMENT HAS BEEN TESTED AND FOUND TO COMPLY WITH THE LIMITS FOR A CLASS A DIGITAL DEVICE, PURSUANT TO PART 15 OF THE FCC RULES.

THESE LIMITS ARE DESIGNED TO PROVIDE REASONABLE PROTECTION AGAINST HARMFUL INTERFERENCE WHEN THE EQUIPMENT IS OPERATED IN A COMMERCIAL ENVIRONMENT. THIS EQUIPMENT GENERATES, USES, AND CAN RADIATE RADIO FREQUENCY ENERGY AND, IF NOT INSTALLED AND USED IN ACCORDANCE WITH THE INSTRUCTION MANUAL, MAY CAUSE HARMFUL INTERFERENCE TO RADIO COMMUNICATIONS. OPERATION OF THIS EQUIPMENT IN A RESIDENTIAL AREA IS LIKELY TO CAUSE HARMFUL INTERFERENCE IN WHICH CASE THE USER WILL BE REQUIRED TO CORRECT THE INTERFERENCE AT HIS OWN EXPENSE.

THIS DIGITAL APPARATUS DOES NOT EXCEED THE CLASS A LIMITS FOR RADIO NOISE EMISSIONS FROM DIGITAL APPARATUS AS SET OUT IN THE INTERFERENCE-CAUSING EQUIPMENT STANDARD ENTITLED "DIGITAL APPARATUS", ICES-003 OF THE DEPARTMENT OF COMMUNICATIONS.

CET APPAREIL NUMERIQUE RESPECTE LES LIMITES DE BRUITS RADIOELECTRIQUES APPLICABLES AUX APPAREILS NUMERIQUES DE CLASSE A PRESCRITES DANS LA NORME SUR LE MATERIEL BROUILLEUR: "APPAREILS NUMERIQUES", NMB-003 EDICTEE PAR LE MINISTRE DES COMMUNICATIONS.

Unauthorized changes or modifications to this equipment are not permitted.

## **TABLE OF CONTENTS**

| INTRODUCTION                         | 4 |
|--------------------------------------|---|
| INSTALLATION                         | 4 |
| USING MENUS TO CONFIGURE THE BALANCE | ŝ |
| PRINT MENU                           | 7 |
| Print Menu Protection                | 7 |
| Reset to Factory Defaults            | 3 |
| Baud Rate                            | 9 |
| Data Bits                            | 9 |
| Parity 10                            | C |
| Stop Bits                            | ) |
| Auto Print Feature1                  | 1 |
| Print Stable Data Only1              | 1 |
| Print Numeric Data Only12            | 2 |
| End 12                               | 2 |
| MENU LOCK-OUT PROTECTION             | 3 |
| USING THE INTERFACE                  | 4 |
| Hardware14                           | 4 |
| RS232 Commands                       | 4 |
| CARE AND MAINTENANCE                 | 7 |
| TROUBLESHOOTING                      | 7 |
| Error Codes1                         | 7 |
| SERVICE INFORMATION                  | 7 |
| I IMITED WARRANTY 19                 | ą |

## INTRODUCTION

The RS232 Interface Accessory is a bidirectional interface which enables either a Precision Standard or an Analytical Standard balance to communicate with a printer or computer equipped with an RS232 serial port. The interface assembly is easily installed in the balance and communication parameters are configured through the balance's front panel.

#### INSTALLATION

Use the following procedure to install the RS232 Interface accessory.

#### **CAUTION**

TO PREVENT DAMAGE TO THE BAL-ANCE, BE CERTAIN THAT POWER IS DISCONNECTED BEFORE REMOVING THE COVER.

1. Disconnect power to the balance.


Remove the draft shield (if in-<sup>1</sup> stalled).

Remove the pan and pan support.

4. Using a philips screwdriver, remove the three screws and washers which secure the cover to the balance. The balance will have to be turned over or on its side to access the screws as they are located on the bottom; two under the front panel and one on the rear.

**CAUTION** 

HOLD THE COVER AND BASE TO-GETHER WHILE REMOVING SCREWS. DO NOT ALLOW COVER OR BASE TO FALL.

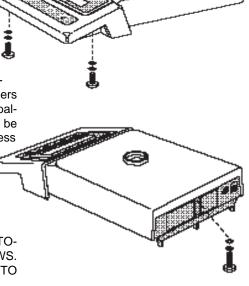



**63** 

- Remove the cover.
- 6. On the rear of the balance, a protective thin plate covers the hole for the 9-pin RS232 connector.

Peel off the plate to expose the hole.

- Place the circuit board and connector assembly in the balance so that the 9pin connector fits through the hole.
- 8. Fasten the assembly using the two screws provided.
- Slide the ribbon connector onto the edge connector of the main circuit board.

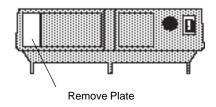


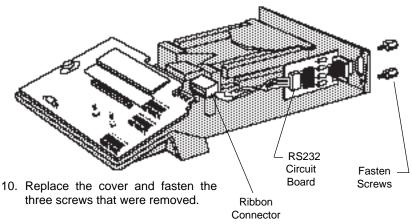

- 11. Replace the draft shield.
- 12. Replace the pan and pan support.
- 13. Reconnect power to the balance.

## **USING MENUS TO CONFIGURE THE BALANCE**

The Analytical Standard and Precision Standard balances both contain four display menus which enable you to calibrate and configure the balance for your specific operating requirements. The print menu shown below is used to configure the RS232 interface.

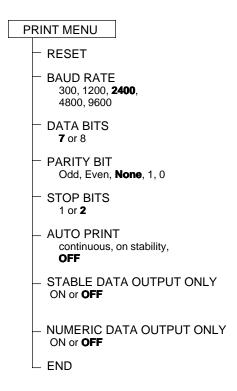

## **PRINT MENU**

When the interface is installed, the Print menu is used to configure the RS232 interface parameters and customize the balance's print functions for your requirements. The following table shows the sequence in which submenus appear on the Print menu.

#### PRINT MENU TABLE

| * r E 5 E Ł | Sets all submenus below to original factory default settings. |
|-------------|---|
| bhud        | Specifies baud rate.  |
| 4AFA        | Specifies number of data bits.                                |
| PAr 1E9     | Specifies parity type, if any.                                |
| 5toP        | Specifies number of stop bits.                                |
| Autop       | Enables/disables Auto print feature.                          |
| SEAPPE      | Enables/disables printing stable-data-only feature.           |
| UU          | Specifies numeric-only or full display data for output.       |
| End         | Used to exit the Print menu and store your selections.        |

Does not appear in menu if menu is locked out.

#### **Print Menu Protection**

The Print menu may be locked out to prevent unauthorized personnel from changing settings. To lock out the print menu, refer to the section titled Menu Lock-Out-Protection.

To access the Print menu press and hold



TARE) until PRINT is displayed, then re-

lease it.

If SAFE is displayed, the Print menu has been locked-out. Settings may be viewed but not changed. See the Menu Lock-Out Protection section to enable it for making changes.

Pr int

To access a submenu:

Repeatedly press MODE until the desired submenu is displayed.

Press TARE to select the displayed submenu.

**NOTE:** You must use END to store any changes you make to the Print menu.

The following sections describe each item on the Print menu in detail.

## **Reset to Factory Defaults**

This submenu enables you to reset all RS232 menu selections to the original factory default settings outlined in the adjacent table.

To reset to factory defaults:

- Access the Reset submenu to view the current setting.
- 2. Press OFF to change the setting.

(Select YES to reset settings or, NO to leave current settings.)

Press (ON to accept the displayed setting.

rESEŁ

| PRINT MENU<br>Factory Defau |         |
|-----------------------------|---------|
| Baud rate                   | br 2400 |
| Data Bits                   | 7 data  |
| Parity                      | None    |
| Stop Bits                   | 2 stop  |
| Auto Print                  | OFF     |
| Stable Data Only            | OFF     |
| Numeric Data Only           | OFF     |

yES no

#### **Baud Rate**

2.

This submenu is used to select the desired baud rate. There are five available baud rates to choose from: 300, 1200, 2400, 4800 and 9600.

bhud

To view or change the baud rate:

- 1. Access the Baud submenu to display the current setting.
- br2400

Press MODE to change the setting. Press (ON to accept the displayed

setting.

ON is released, BAUD will be When ( displayed again and the Print menu will be returned.

# **JAFA**

#### **Data Bits**

The total number of bits for Data, Parity and Stop must equal 9 or 10. (see examples). The balance will not permit you to select a combination that does not equal 9 or 10.

To set the number of data bits to 7 or 8:

- Access the Data submenu to display the current setting.
- Press (MODE) to change the setting. 2.

Press (ON to accept the displayed setting. When (on take) is released, DATA will be displayed again and the Print menu will be returned.

#### **EXAMPLES**

8 Data + 2 Stop + No Parity

8 Data + 1 Stop + Odd Parity = 10

7 Data + 1 Stop + Odd Parity = 9

7 dafa

8 4868

## **Parity**

Parity can be set to Odd, Even, None, or a marker of 0 or 1 as follows:

- 1. Access the Parity submenu to display the current setting.
- 2. Press (MODE) to change the setting.

Press to accept the displayed setting.

**NOTE:** If all selections do not appear, total number of data, parity and stop bits is currently < 8 or > 10. Data or stop bits must be changed.

When TARE is released, PARITY will be displayed again and the Print menu will be returned.

## Stop Bits

The number of stop bits can be set to 1 or 2 as follows:

- 1. Access the Stop submenu to display the current setting.
- 2. Press (MODE) to change the setting.
- 3. Press (ON to accept the displayed setting.

**NOTE:** If all selections do not appear, total number of data, parity and stop bits is currently < 8 or > 10. Data or parity bits must be changed.

When TARE is released, STOP will be displayed again and the Print menu will be returned.

PAr ity

nanE

-044-

- 1-

-8-

SEOP

I SEOP

2 SEOP

#### **Auto Print Feature**

When enabled, the Auto Print feature causes the balance to automatically output display data in one of two ways: continuously, or upon stability.

AutoP

To select one of these Auto Print methods, or to turn the feature off:

Access the AutoP submenu to display the current setting.

OFF

2. Press NODE to change the setting.

Cant

Press (ON to accept the displayed setting.

On 566

When on is released, AUTOP will be displayed again and the Print menu will be returned.

## **Print Stable Data Only**

When enabled, this feature permits only stable display data to be output. To set the feature ON or OFF:

SEABLE

- Access the Stable submenu to display the current status.
- 2. Press OFF to change the status.

On

Press (N) to accept the displayed status.

OFF

When TARE is released, STABLE will be displayed again and the Print menu will be returned.

## **Print Numeric Data Only**

This submenu is used to select numeric data only, or full display data for RS232 output. Set this feature ON to output numeric display data only, or OFF to output full display data as follows:

UR

- 1. Access the Nu submenu to display the current status.
- 2. Press MODE to change the status.

Press (TARE) to accept the displayed status.

0n

When tare is released, NU will be displayed again and the Print menu will be returned.

OFF

#### **End**

You must use END to exit the Print menu. Changes you make in the Print menu are only stored in memory if you use End.


End

To exit the Print menu and store your selections, press (NTARE) when END is displayed. The balance will be returned to normal weighing operations.

## **MENU LOCK-OUT PROTECTION**

Access to the Print Menu can be disabled using the lock-out switch located under the right side of the balance, near the display.

- 1. Turn display off and unplug the power cord.
- Slide the balance toward you, with the front over the edge of a table. (You can also turn the balance on its left side, but if you do, you MUST remove the draft shield, pan and pan support first!)
- Locate hole under display where switch is located.
- Using a small screwdriver, slide the switch forward for LOCKED or back for UNI OCKED.
- 5. Plug in the power cord and turn on the balance.

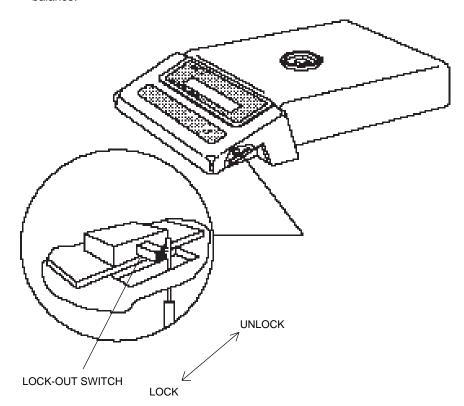

## USING THE INTERFACE

When the interface is installed, a short



press on the MODE switch will initiate a

print command. Data output is in the format shown under the P command in the RS232 Command Table.

When the interface is connected to a computer, two way communication between the computer and balance is possible using the commands outlined in the RS232 Command Table.

#### **Hardware**


The balance can be interfaced with other equipment utilizing the 9-pin subminiature "D" connector. The pinout and pin connections are shown in the adjacent illustration.

The balance will not output any data unless pin 5 (CTS) is held in an ON state (+3 to +15 VDC). Interfaces not utilizing the CTS handshake may tie pin 5 to pin 6 to defeat it.

#### **RS232 Commands**

All communication is accomplished using standard ASCII format. Only the characters shown in the following table are acknowledged by the balance. Any other commands, control characters or spaces are ignored.

Commands sent to the balance must be terminated with a carriage return (CR) or carriage return-line line feed (CRLF). For example, a tare command should appear as shown in the adjacent diagram. Data output by the balance is always terminated with a carriage return - line feed (CRLF).

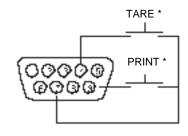

- 5VDC (50 mA max.) 1
- 2 Data Out (TXD)
- 3 Data In (RXD)
- Tare (External signal)
- Clear To Send (CTS) 5
- 6 Data Terminal Ready (DTR)
- Ground 7
- Request To Send (RTS)
- Print (External signal)
- External PRINT and/or TARE switches may be installed as shown in the diagram. Momentary contact switches must be used.

#### TARE COMMAND

Field: Length:

| Т | CR | LF |
|---|----|----|
| 1 | 1  | 1  |

## **RS232 COMMAND TABLE**

| Command<br>Character | Description  |   |
|----------------------|--|---|
| ?                    | Print current mode   | Field: Mode Stab CR LF Length: 5 1 1 1  blank if stable "?" if unstable  Grams Momme Pennyweight Pounds Carats Pounds:ounces Avoidupois ounces Custom unit Troy ounces Parts counting Grains Percent weighing                 |
| nnnA                 | Set Auto Print feature to "n" (see table).   | Taels Error  n = 0 Turns feature OFF n = S Output on stability n = C Output is continuous   |
| С                    | Begin span calibration   |   |
| хD                   | Set 1 second print delay (set x  | x = 0 for OFF, or $x = 1$ for ON)   |
| ХĬ                   | Set Averaging Level to "x", where $x = 0$ to 3 (see table).  | 0 = minimum level<br>1 =<br>2 =<br>3 = maximum level  |
| L                    | Begin linearity calibration  |   |
| P                    | Print display data  When "numeric only" display data is selected for output in the RS232 menu, the Mode field is not output. | Field: Weight Length: 9 1 5 1 1 1 1  Same as? command Displayed weight sent right justified w/lead zero blanking.  Nine characters include: decimal point (1) weight (7 max)) polarity (1): blank if positive "-" if negative |

## **RS232 COMMAND TABLE (Cont.)**

| Command<br>Character | Description   |  |  |
|----------------------|---|--|--|
| xS                   | Set stable data only printing (set $x = 0$ for OFF, or $x = 1$ for ON).                         |  |  |
| Т                    | Same effect as pressing tare button   |  |  |
| V                    | Print EPROM version  Field: Model # EPROM # CR LF 7 15 1 1  Balance Model  "98101-XX Sr*XX.X"   |  |  |
| xΖ                   | Set Auto Zero to "x", where x = 0 to 3 (see table).  0 = OFF 1 = .5 d 2 = 1 d 3 = 3 d           |  |  |
| Esc L                | Prints listing of Setup and Print menu settings.  |  |  |
| Esc R                | Resets Setup and Print menus to factory defaults. CAUTION: This will reset RS232 configuration. |  |  |
| Esc S                | Save current settings.  |  |  |

## **CARE AND MAINTENANCE**

The RS232 Interface Accessory once installed does not require maintenance.

## **TROUBLESHOOTING**

| SYMPTOM                      | PROBABLE CAUSE(S)                        | REMEDY  |
|------------------------------|--|---|
| RS232 interface not working. | Print menu settings not properly set up. | Verify interface settings in Print menu correspond to those of peripheral device. |
|                              | Cable connections.                       | Check cable connections.  |

#### **Error Codes**

When the RS232 interface is installed in Precision Standard and Analytical Standard balances, the following error codes are added to the standard list of error codes:

- 4.0 Bad RS232 frame. Check RS232 menu parameters and correct.
- 4.4 RS232 buffer is full. May occur if no printer or computer is connected to the interface. To clear buffer turn balance off, or enter and exit Print menu and select FND.
- 4.5 Function is disabled by the Lockswitch.

## **SERVICE INFORMATION**

If the Troubleshooting section does not resolve or describe your problem, you will need to contact an authorized Ohaus Service Agent. For Service assistance in the United States, please call Ohaus Corporation toll-free at (800) 526-0659. An Ohaus Product Service Specialist will be available to help you.

#### LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.



Ohaus Corporation 29 Hanover Road, Florham Park, NJ 07932, USA Tel: (201) 377-9000,

Fax: (201) 593-0359

Ohaus Europe Ltd. ENGLAND Tel: +44 (0) 1954 251343, Fax: +44 (0) 1954 250205

With offices in Germany, France, Spain, Italy, Canada, Mexico and Japan.

P/N 77017-01 R0296 © Ohaus Corporation 1993, 1996 all rights reserved.